


Installation Instructions: MATLAB Cache Re-Location & BAX® Real-Time *Salmonella* Patch


Software Version Verification

1. Verify the BAX® System Q7 Software version is 3.6.7262
 - a. Go to the “Help” menu
 - b. Select the “About BAX® System” option
 - c. Once version is verified, continue to download and install instructions
 - i. If version is different from above, contact Diagnostics Support, diagnostics.support@hygiena.com for assistance or upgrade to the latest software version located [here](#)

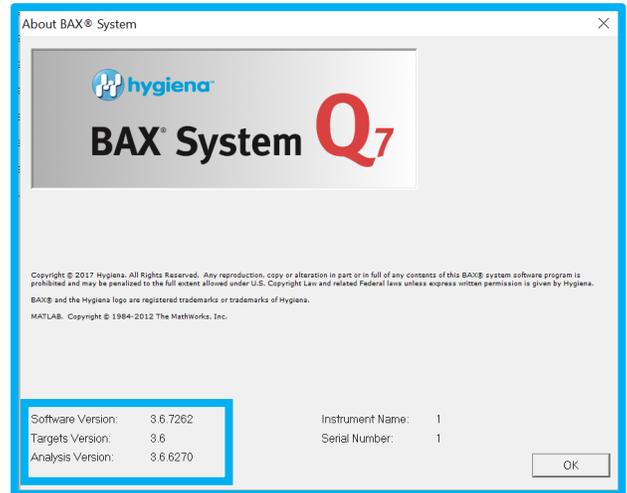


MATLAB Cache + BAX® Real-Time *Salmonella* Patch

1. Using the link below, download the **3.6.7032.1383.zip** folder to a known location
 - a. <https://www.hygiena.com/bax-system/q7-software-improvement.html>
 - b. Optional: Download to a memory stick/flash drive to transport zip file to BAX® System Q7 computer
2. Unzip the downloaded .zip file in the known location
 - a. Right-click the .zip file and select “Extract All”
 - b. Unzip the file to any known location (e.g. downloads, desktop, etc.)
3. Open the uncompressed folder and run the **Update.bat program as an administrator**
 - a. Right-Click the Update file and select “Run as Administrator” (if this option is not available you will need to contact your IT department to install the patch)
 - b. Follow the command prompts
 - c. Analysis version is installed after finishing prompts
 - d. Hit any key to close the Command window.
 - e. Re-check the software version to ensure proper installation by following section 1 above.

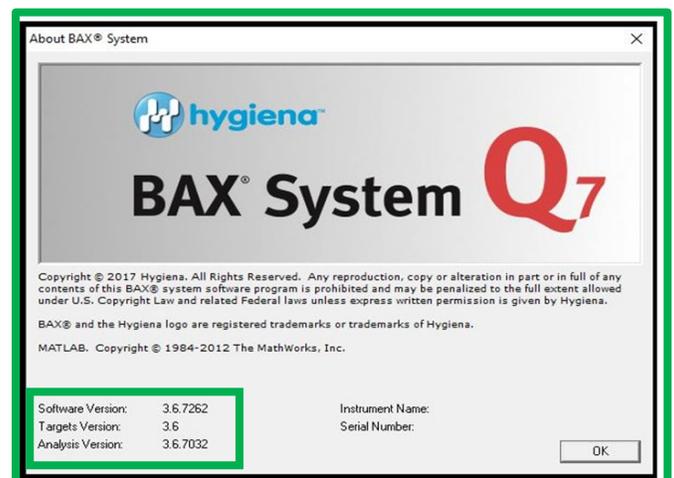

MATLAB Cache ONLY

1. Using the link below, download the **CacheMove.zip** folder to a known location
 - a. <https://www.hygiena.com/bax-system/q7-software-improvement.html>
 - b. Optional: Download to a memory stick/flash drive to transport zip file to BAX® System Q7 computer
2. Unzip the downloaded .zip file in the known location
 - a. Right-click the .zip file and select “Extract All”
 - b. Unzip the file to any known location (e.g. downloads, desktop, etc.)
3. Open the uncompressed folder and run the **Update.bat program as an administrator**
 - a. Right-Click the Update file and select “Run as Administrator” (if this option is not available you will need to contact your IT department to install the patch)
 - b. Follow the command prompts
 - c. Analysis version is installed after finishing prompts
 - d. Hit any key to close the Command window.
 - e. Re-check the software version to ensure the analysis version did not change since only the MATLAB cache relocation script was installed

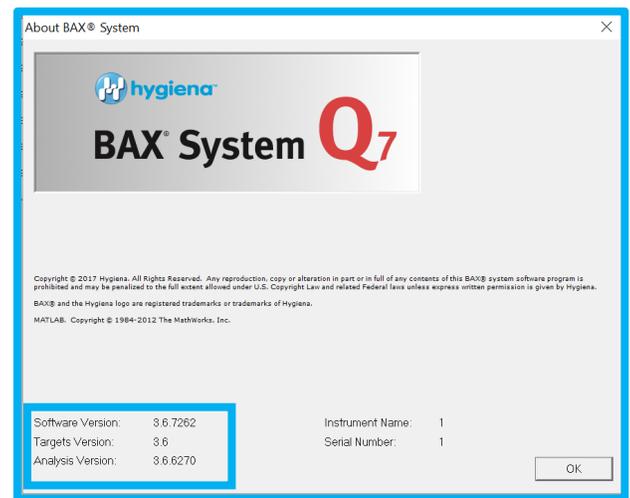

Software Revert

1. If there is a need to revert to the previous software, utilize the **Revert.bat or DeleteMCREnv.bat** program located in the same folder.
2. After running the reversion program, you must **restart the computer**.
3. After restart, re-check the software version by following the instructions in the software version verification section.